

Millennia Pro Systems Pre-Installation Guide



Spectra-Physics, Inc.
3635 Peterson Way
Santa Clara, CA 95054
April 2020

Table of Contents

Introduction.....	3
When the system Arrives	4
Inspection	4
Review Instruction Manuals.....	4
Laser Safety Considerations.....	4
Diagnostics	4
Pre-Installation Considerations	5
Location and Environment.....	5
Physical Description	6
Dimensions and Weights.....	6
Utility Requirements	6
Power Requirements.....	6
Pre-Installation Checklist	7
Physical Locations:.....	7
When Your System Arrives:.....	7
Maintenance Agreement.....	8
Nalco Mixture	8
Filters.....	8
Maintenance Agreement (Continued)	9
Chiller Fluid	9
General Procedures	9

Introduction

Congratulations on your purchase of a Spectra-Physics system. This guide describes pre- installation information for your laser system.

Some preliminary planning is essential to avoid unnecessary delays during installation and to ensure easy operation and access to your system. You are requested to carefully consider your operating environment prior to installation. Proper water, power and room temperature are required for each system.

A checklist of pre-installation considerations is provided in this guide. You are responsible for meeting these requirements prior to installation, with due consideration given to all applicable building and safety codes.

We at Spectra-Physics intend to provide you with responsive support so that you can derive great satisfaction and value in using our systems for your applications. We are available to you at 1-800-456-2552.

When the system Arrives

Inspection

When the system arrives, any sign of damage to the shipping crates should be brought to the attention of the delivering freight company. A claim must be filed with that commercial carrier (usually within 30 days). Notify the originating Spectra-Physics office of any shipping damage. Shipping damage is not covered by Spectra-Physics.

Your packing list will show all items that you have ordered. Open all the packages and check each item for possible damage during shipping. Check the items against your packing list. Some items may have been installed at the factory.

Each system comes with a manual; verify that you have received this item.

Please report any missing or damaged items to Spectra-Physics, or you may contact your Spectra-Physics Sales Engineer.

Review Instruction Manuals

Please read the manual to get vital information about your system. Familiarize yourself with the system. You are encouraged to spend as much time as possible reviewing the system components before your Spectra-Physics Service Engineer arrives for the installation and training.

Laser Safety Considerations

In addition to reviewing the sections in the manual regarding laser safety; be sure to have the proper safety glasses available for ALL lab personnel present during the installation and testing of your system. For more information, please call Spectra-Physics at 1-800-456-2552.

Diagnostics

During the course of installation, power measurements will be demonstrated on all of the appropriate wavelengths. Your Service Engineer can identify which specifications will be demonstrated and the equipment necessary to conduct such tests. A non-standard system will require special consideration.

To have other published specifications demonstrated, consult with your Service Engineer to determine whether additional diagnostic equipment will be required.

Pre-Installation Considerations

Location and Environment

The location of the system and environment of your lab should have the following features:

- A. A safe location that meets all applicable building codes.
- B. Easy access with adequate clearance around the instrument.
- C. An optical table that will meet the space requirement of the instruments to be installed.
- D. Proper air conditioning could be critical for the performance of the laser. For certain applications ambient room temperature changes may be an important factor for the laser system's performance. Air ducts should not blow directly on laser or optical path.
- E. To ensure stable day-to-day operation, the recommended minimum and maximum operating room temperatures are 18 - 35°C. Room temperature should ideally be 22°C. See additional specifications in the User's Manual.
- F. In some applications vibration isolation may be required for your system. Structural integrity of the flooring could play an important role.
- G. Please note that the laser head requires a sturdy and stable optical table.

Consider room requirements for future maintenance and upgrades by your Spectra- Physics Field Service Engineer.

Physical Description

Dimensions and Weights

Feature	Specifications	
Size (W x H x D)	English Unit	Metric Unit
Laser head	5.3 x 4.1 x 14.0 in.	13.5 x 10.4 x 35.0 cm
Power supply	19.0 x 6.9 x 17.9 in.	48.3 x 17.5 x 45.5 cm
Umbilical length	10 ft.	3 m
Weight		
Laser head, scientific	22.0 lb.	10 kg
Power Supply (J40)	55.0 lb.	25.0 kg
Power Supply (J80)	58.0 lb.	26.4 kg
Thermoroack Chiller	39.0 lb.	18 kg

Utility Requirements

Water

The Millennia system is a closed loop water cooled system and requires no utility water services. See User Manual for additional information.

Fill the chiller reservoir with distilled water. Do not use deionized water and do not place the chiller above the laser. Should the unit not be installed properly, and a leak develops, dripping water may damage the laser. The chiller must always be on when the power supply is on, even if the diode lasers are not on.

Power Requirements

Electrical

Power Supply Single Phase¹
 100 to 240 Vic, 50/60 Hz, T 6.3 A, 250 V fuses (two)

Thermorack chiller

Universal Input 100-240 VAC, 50-60 Hz, 9-4 amps max

Two modes of control are offered for the standard Millennia system:

- A. A handheld Controller provides local control of the system via an 8-foot cable that connects to the front of power supply. For more information on the controller, refer to the User's Manual.
- B. For users who prefer to operate the laser remotely, either directly or via a computer program, a standard RS232 serial link is provided on the power supply to connect to a computer or terminal.

Both methods are described in Chapter 6 of the User's manual.

Pre-Installation Checklist

Before the arrival of your Spectra-Physics' Service Engineer please review the following pre- installation requirements. When all the requirements have been met, initial the boxes and fax a signed copy to the Spectra-Physics Service department at **(408) 980-6921**.

Physical Locations:

- A location with adequate clearance around system to conduct service and accessible by Spectra-Physics personnel.
- A temperature-controlled room.
- Utility services have been installed.
- Local building and safety codes are in compliance and have been verified.

When Your System Arrives:

- Check crates for damage.
(If damaged, file a claim with the carrier and notify Spectra-Physics.)
- Uncrate and place the system on your work surface.
(Two or more people may be required to lift some equipment.)
- Compare the packing list with your quotation. Call your Spectra-Physics office about any discrepancies.
- Check that all manuals were received.
- Save all packing and shipping material until the installation has been completed.
- Obtain the correct safety glasses and a power meter.

Customer Signature

Date

Phone Number

Fax Number

E-Mail

Sales Order Number

Maintenance Agreement

To maintain a valid warranty on your Spectra-Physics system, it is necessary for the customer to assume the responsibility and perform the routine maintenance program. Failure to do so may result in the warranty being voided.

Nalco Mixture

The Nalco Cleaner (P/N 1607-0547) is used to flush your system clean of ferrous metals and copper alloys as a result of corrosion. It is recommended as a maintenance procedure that the chiller/critical cooling loop in each piece of equipment is cleaned and treated 1x per year. The laser does not have to be lasing. The cleaner should be circulated for a minimum of 8 hours (the longer the better if time permits in order to assure thorough cleaning).

The Nalco Solution (P/N 1607-0546) is a premixed liquid corrosion inhibitor designed for use in closed loop cooling systems. It can be added directly to a closed cooling system (used in place of water); resulting in a proper treatment without dilution for up to 6 months (even if the pink tracer turns clear). This solution is a complete inhibitor that protects ferrous metals and copper alloys from corrosion. It is nitrite free and minimizes the challenge of bacteria control (depending on environment and usage).

Filters

There is 10 Micron Particle Water Filter located at the rear of the chiller that should be inspected every 3 months and changed when dirty or no longer able to rinse clean. Refer to User manual for additional info. You should also monitor the laser tower temperature to make sure cooling is within specifications.

Water filter replacements can be ordered using P/N 2604-0490.

There is also an Air filter that is located behind the right-hand grill on the power supply front panel (J40 only). Refer to User Manual for additional info. It should also be inspected and blown clean every 3 months.

Air filter replacements can be ordered using PN 0129-9322S.

If the laser is not going to be used for an extended period of time, turn off the laser and drain the water or cooling solution.

Regularly inspect filter on the power supply for any buildup of dust. If dirty, remove the filter and wash it in clean water. The rate at which the filter becomes more restrictive is mainly dependent on the cleanliness of the environment.

Maintenance Agreement (Continued)

***NOTE:** to restrict the growth of algae in the reservoir, it is recommended that the reservoir cover be kept in place and that all circulation lines be opaque. This will eliminate the entrance of light that is required for the growth of most common algae.

Chiller Fluid

The chiller fluid level should be appropriate for the model laser and must meet Spectra-Physics specification requirements. Insufficient cooling could result in a decrease of performance or damage to the laser system.

General Procedures

The chiller operation of the laser system should be checked once a week. The power output of the system should also be checked once a week by verifying the output values.

If any of the output characteristics have changed, please call Spectra-Physics Technical Support at **1-800-456-2552**.