# Mai Tai<sup>®</sup>

## Ultrafast One Box Ti:Sapphire Lasers

# • MKS | Spectra-Physics

The Spectra-Physics Mai Tai Ti:Sapphire oscillator is used for a variety of applications ranging from general spectroscopy to multiphoton microscopy. With its leading performance specifications and largest installed base of any laser of its type, the Mai Tai is truly a cutting-edge tool for basic research and biological imaging.

The Mai Tai series includes three models of increasing power and tuning range to allow customers to choose the system to best match their application. Ideal for multiphoton microscopy, the Mai Tai HP provides more than 300 nm (690–1040 nm) in useable tuning range with over 2.5 W of average power and a pulse width of less than 100 fs. For customers interested in systems with a shorter pulse width, the Mai Tai BB provides more than 240 nm in useable tuning range with a pulse width of less than 80 fs. Finally, the Mai Tai XF-1 provides more than 200 nm in useable tuning range with an even shorter pulse width of less than 70 fs. For access to longer wavelengths beyond 1040 nm, Spectra-Physics offers InSight<sup>®</sup> X3<sup>™</sup> with exceptionally wide tuning range (680–1300 nm) for deep in vivo imaging.

The Mai Tai series reliability is maintained using proven, ultra-stable regenerative mode-locking. Using this method, the Mai Tai oscillator is capable of hands free, drop-out free wavelength tuning enabling speedy collection of excitation profiles all at the click of a mouse. StabiLok<sup>®</sup> real-time monitoring technology also enhances system reliability by providing excellent beam pointing stability and minimal average power fluctuations, as well as eliminating wavelength drift. These two features combine to make the Mai Tai oscillator the most reliable and most versatile hands-free laser source available.

### The Mai Tai Advantage

- Short output pulse width to provide high peak power
- Wide tuning range for efficient excitation of all commonly used fluorophores
- Smooth tuning with no pulse drop outs
- Outstanding pointing stability eliminates the need to realign experimental set-up
- Regenerative mode locking for unmatched stability

#### **Applications**

- Multiphoton microscopy
- Time-resolved photoluminescence
- Non-linear spectroscopy
- Optical computed tomography
- Surface second harmonic generation
- Terahertz imaging
- Semiconductor metrology
- Materials processing
- Amplifier seeding

# Mai Tai Specifications<sup>1, 10, 11</sup>

	Mai Tai HP	Mai Tai BB	Mai Tai XF-1					
Output Characteristics								
Pulse Width <sup>2, 3</sup>	<100 fs	<80 fs	<70 fs					
Tuning Range <sup>4</sup>	690–1040 nm	710–990 nm	710–920 nm					
Average Power at 800 nm <sup>2</sup>	>2.5 W	>1.5 W	>900 mW					
Average Power, Alternative Wavelengths⁵	>500 mW at 690 nm >1.35 W at 710 nm >1.35 W at 920 nm >300 mW at 1040 nm	>650 mW at 710 nm >650 mW at 920 nm >250 mW at 990 nm	>400 mW at 710 nm >400 mW at 920 nm					
Beam Roundness <sup>2</sup>	0.9–1.1							
Astigmatism <sup>2</sup>	<10%							
Repetition Rate <sup>2, 6</sup>	80 MHz ±1 MHz							
Beam Pointing Stability	<50 μrad/100 nm							
Noise <sup>2, 7</sup>	<0.15%							
Stability <sup>8</sup>	<±1%							
Spatial Mode <sup>2</sup>	TEM <sub>00</sub> , M <sup>2</sup> <1.1							
Polarization	>500:1 Horizontal							
Beam Divergence <sup>2</sup>	<1.2 mrad							
Beam Diameter (1/e2)2	<1.2 mm							
Physical Characteristics								
Size (L x W x H)	23.89 x 13.79 x 5.95 in (606.81 x 350.27 x 151.13 mm)							
Weight – Laser Head	70 lbs (32 kg)							
Umbilical Length	10 ft (3 m)							
Environmental Requirements								
Altitude	Up to 2000 m							
Temperature, Operating	20–25°C							
Relative Humidity, Operating	Maximum 75% non-condensing up to 25°C							
Temperature, Storage	15–35°C							
Relative Humidity, Storage	<65% for 15–35°C							
Cooled Water Temperature in Closed-loop Chiller	21°C typical <sup>9</sup>							

2. Specification applies to 800 nm only.

3. A sech2 pulse shape is used to determine the pulse width as measured with a Newport PulseScout® autocorrelator.

4. Mai Tai is also available with a fixed, factory preset wavelength within the wavelength range noted.

5. Specifications apply to operation at the wavelength noted.

6. Laser operation is specified at a nominal repetition rate of 80 MHz.

7. Specification represents rms noise measured in a 10 Hz to 10 MHz bandwidth.

8. Percent power drift in any 2-hour period with <±1°C temperature change after a 1-hour warm up.

9. Avoid obstructing the air exhaust grills which will result in the recirculation of hot exhaust air. Cooling air enters through the front panel and exits through the rear fan apertures.

10. Mai Tai eHP is available with short pulses.

11. The Mai Tai is a Class IV – High-Power Laser, whose beam is, by definition, a safety and fire hazard. Take precautions to prevent exposure to direct and reflected beams. Diffuse as well as specular reflections can cause severe skin or eye damage.

# • MKS | Spectra-Physics



## Mai Tai Dimensional Drawing





www.spectra-physics.com

MaiTai\_05/25 ©2025 MKS Inc. Specifications are subject to change without notice.

#### 1565 Barber Lane, Milpitas, CA 95035 USA PHONE: 1-800-775-5273 1-408-980-4300 FAX: 1-408-980-6921 EMAIL: sales@spectra-physics.com

Belgium	+32-(0)0800-11 257	belgium@newport.com	Korea	+82-31-8021-1600	korea@spectra-physics.com
China	+86 510 8113 2999	spectra-physics-china@mksinst.com	Netherlands	+31-(0)30 6592111	netherlands@newport.com
France	+33-(0)1-60-91-68-68	france@newport.com	Singapore	+65-6664-0040	sales.sg@newport.com
Germany	/ Austria / Switzerland		Taiwan	+886-3-575-3040	sales@newport.com.tw
	+49-(0)6151-708-0	germany@newport.com	United Kingdom	+44-1235-432-710	uk@newport.com
Japan	+81-3-3556-2705	spectra-physics.ip@mksinst.com			

MKS products may be subject to export, re-export, and economic sanctions controls administered by multiple global jurisdictions and may include the United States. Export, re-export, diversion, transfer, or use contrary to all applicable laws is prohibited. Spectra-Physics<sup>®</sup>, InSight<sup>®</sup>, Mai Tai<sup>®</sup>, StabiLok<sup>®</sup> and PulseScout<sup>®</sup> are registered trademarks, and X3<sup>™</sup> is a trademark of MKS Instruments, Inc. or a subsidiary of MKS Instruments, Inc. Spectra-Physics Milpitas, California, Stahnsdorf, Germany, Rankweil, Austria and Rehovot, Israel have all been certified compliant with ISO 9001.